



# How to fit your eyepiece with a reticle?

**Important notice:** This manual will only show you how to put a reticle in your eyepiece in a Nikon or Olympus microscope. For other microscopes, please contact your local service agent, they can tell you exactly which size reticle you need. During your yearly service check, the technician can fit the reticle in the eye piece for you.



**Step 1:** Take your eye-piece out of the tube and unscrew the small ring at the lower end as indicated by the red eclipsis..

**Step 2:** Now measure precisely the diameter of the eyepiece. There is a small rest where the reticle will be placed on.

In this example, the diameter is 21,4 mm.

The thickness of the reticle for a Nikon microscope will be 1,5 mm and for an Olympus 1,0 mm.



**Step 3:** Carefully place the supplied reticle in the eyepiece and replace the ring taken out during step 1.

Now you will have to calibrate your microscope. A manual is available on [www.leja.nl](http://www.leja.nl)